AK Eychmüller GroupSeminar

Brief introduction to the
WinXPow for XRD.

AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

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TU Dresden » Faculty of Science » Department of Chemistry and Food Chemistry » PC2

GROUP

- Start
- News
- Events
- Members
- WG Bund
- Jobs Vacancies
- Internal Info

RESEARCH

- Nanocrystals
- Projects
- Publications
- Cooperations

TEACHING

- Overview
- Lectures
- Exercises
- Practical Training
- Miscellanea

@akeychm excursion 2007

important links

Information system on hazardous substances

- [Databases on hazardous substances of BGIA](#)
- [material safety data sheet of euSDB](#)

important contact

- [Glassblower](#)

Analysis and Service (Department of chemistry)

- [Overview of measurements and equipment](#)
- [Measuring data service](#)
- [nanoeva](#)

downloads

Seminar - on wednesdays at 9.30 a.m./room 302/E.-Müller-Bau

- [current seminar plan](#)

Ordering and application form

- [Order form of XRD-measurements](#)
- [Chemicals ordering](#)
- [POSTERPRINT-instruction](#)

Safety and Disposal

- [Eye and lab shower controle](#)
- [Solvent Disposal and Refill](#)

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Mailprograms destroy .raw-files!! (Thunderbird, Outlook, etc.)


⇒ downloading from:

1. Measuring Data Service
2. Squirrel Mail

AK Eychmüller GroupSeminar


Brief introduction to the WinXPow for XRD.

- Licenced to:
Professorship of the Inorganic Chemistry
Prof. S. Kaskel
- Permission for us:
installation on one computer for the hole group
⇒ office R 003
⇒ user: office03
⇒ psw: useropen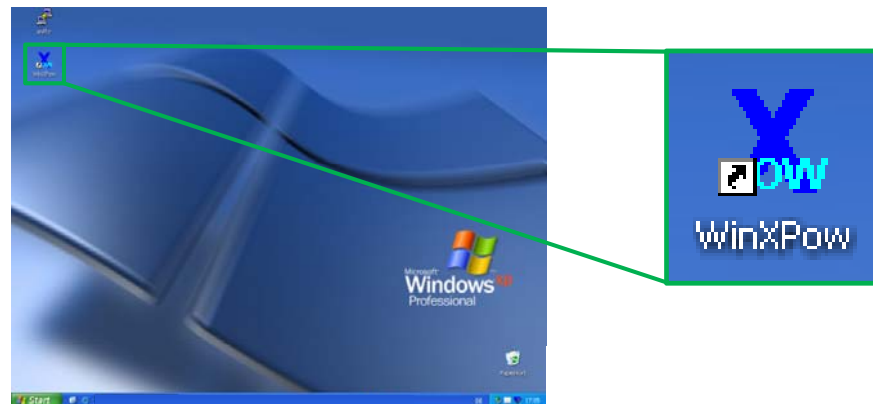

AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

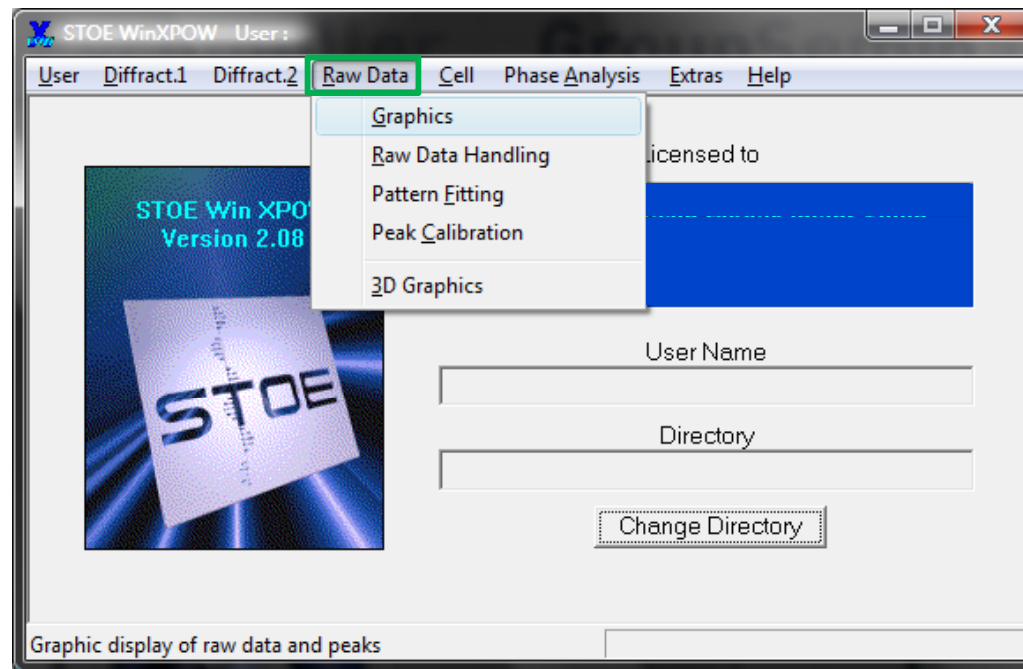
AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Raw Data → Graphics


- File → Open → cw_Alloy-X1.raw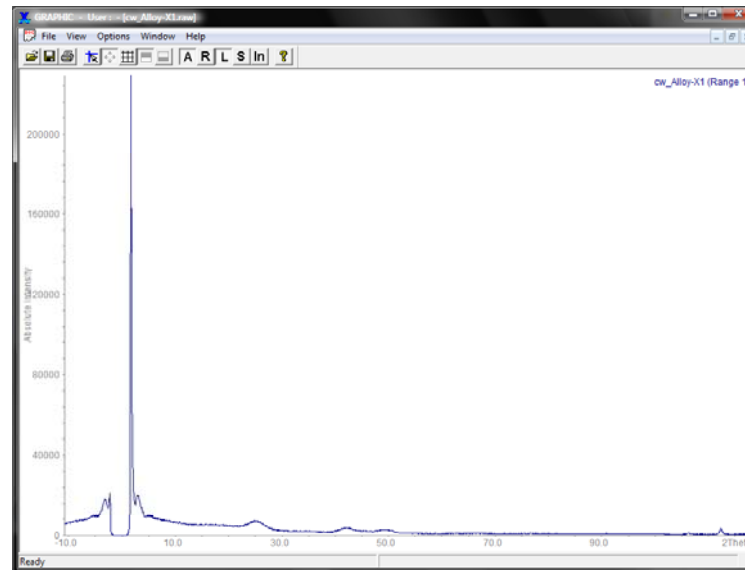

AK Eychmüller GroupSeminar


Brief introduction to the WinXPow for XRD.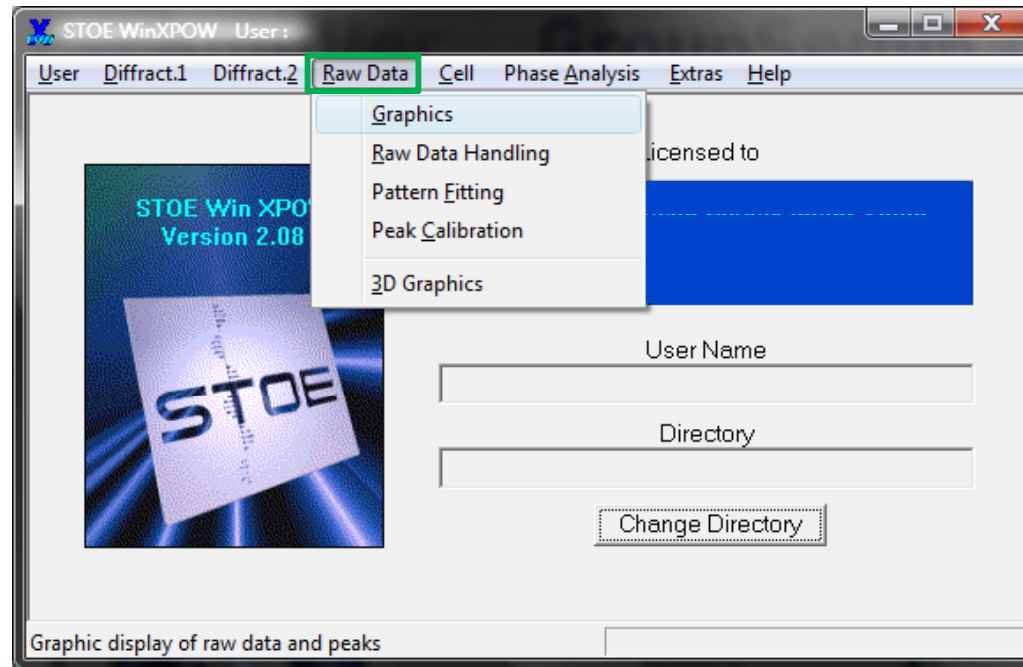


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Raw Data → Raw Data Handling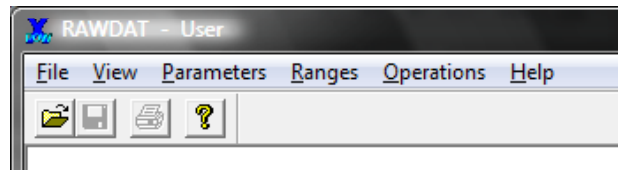


- File → Open → cw_Alloy-X1.raw

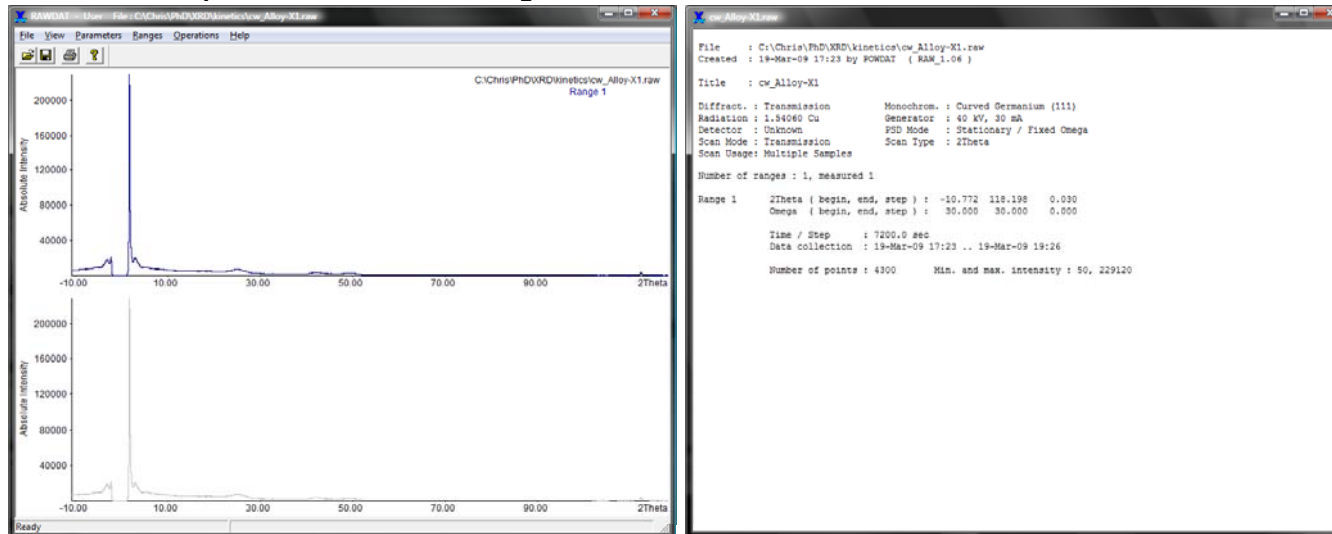AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.


- Ranges → Truncate Range → 2Theta: 10 to 90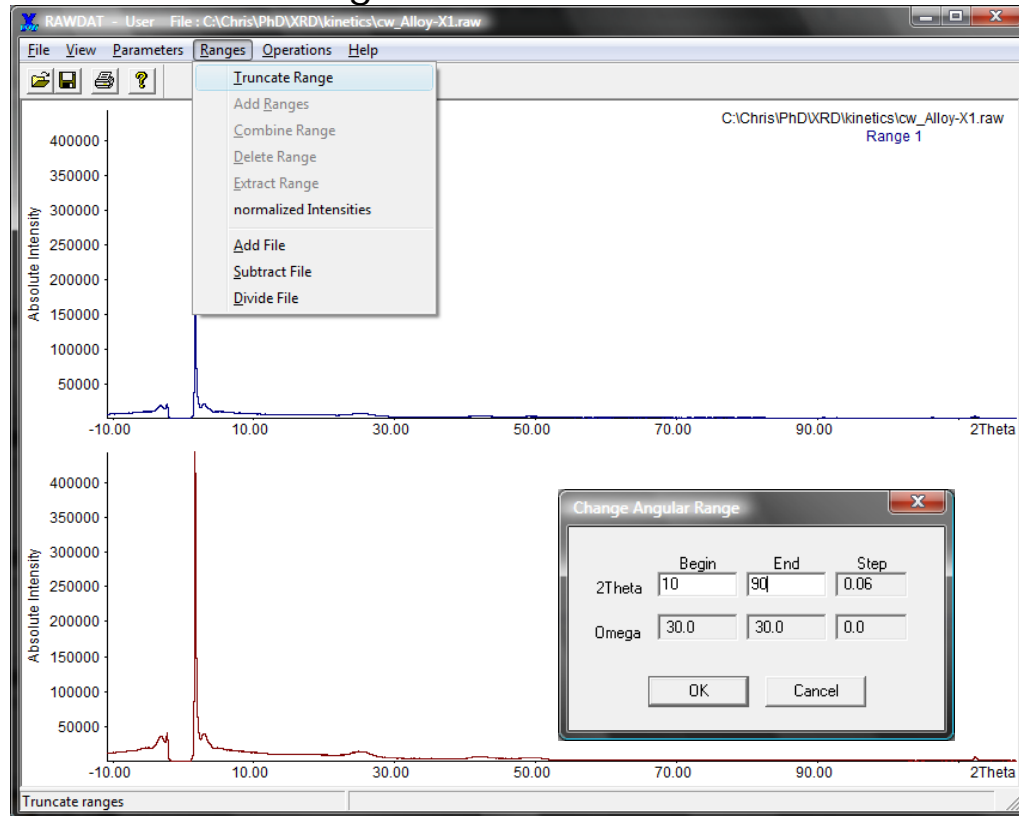

AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.


- File → Save As → cw_Alloy-X1opt.raw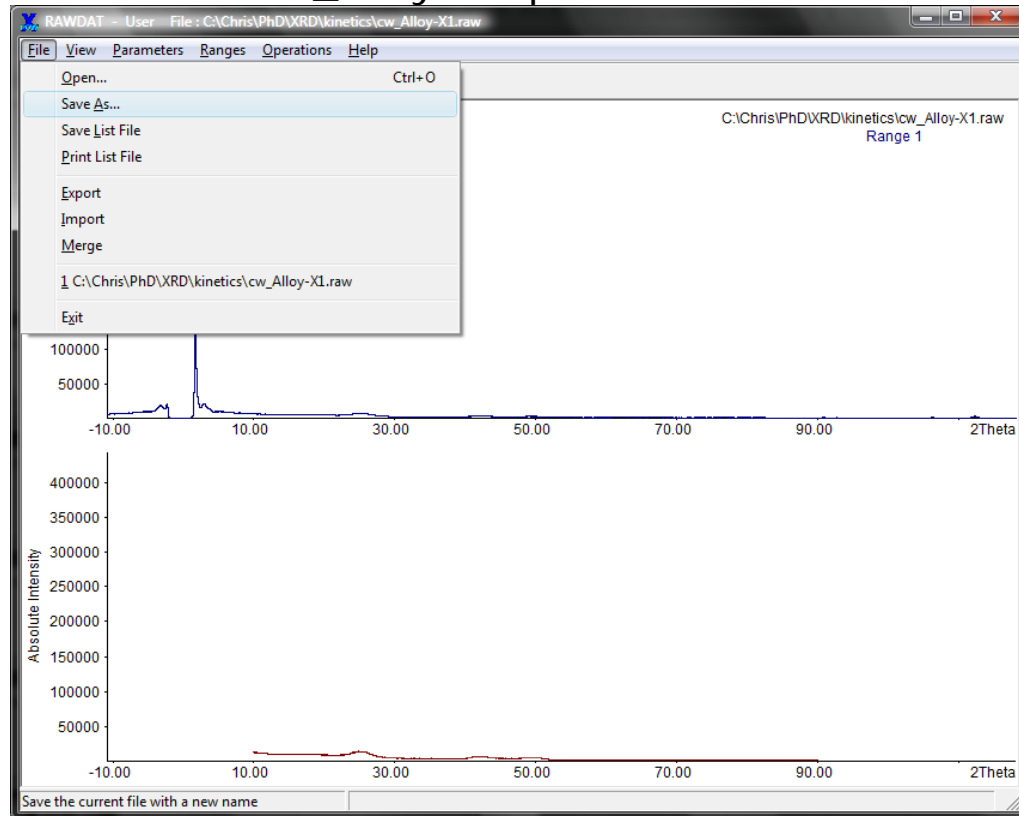


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Raw Data → Graphics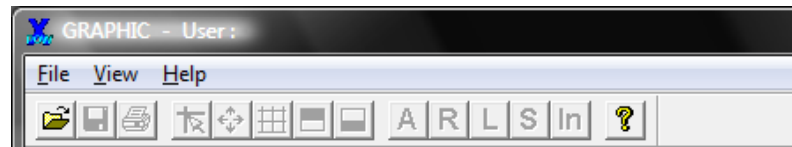


- File → Open → cw_Alloy-X1opt.raw

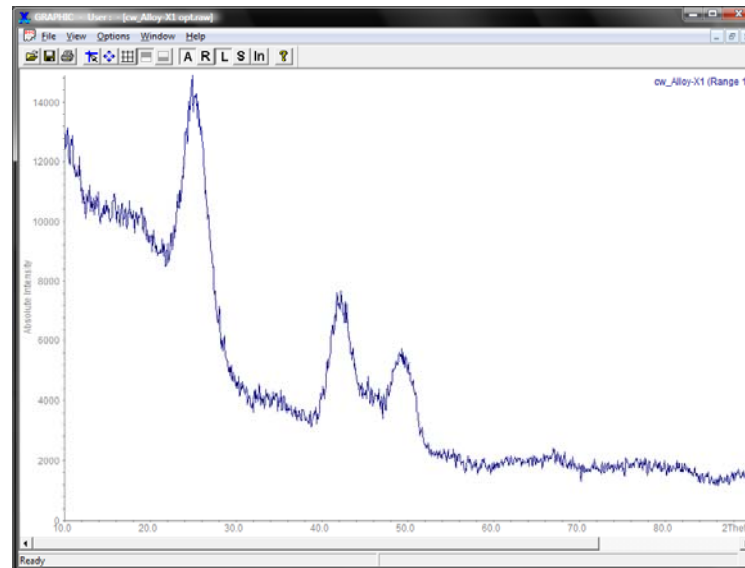AK Eychmüller GroupSeminar


Brief introduction to the WinXPow for XRD.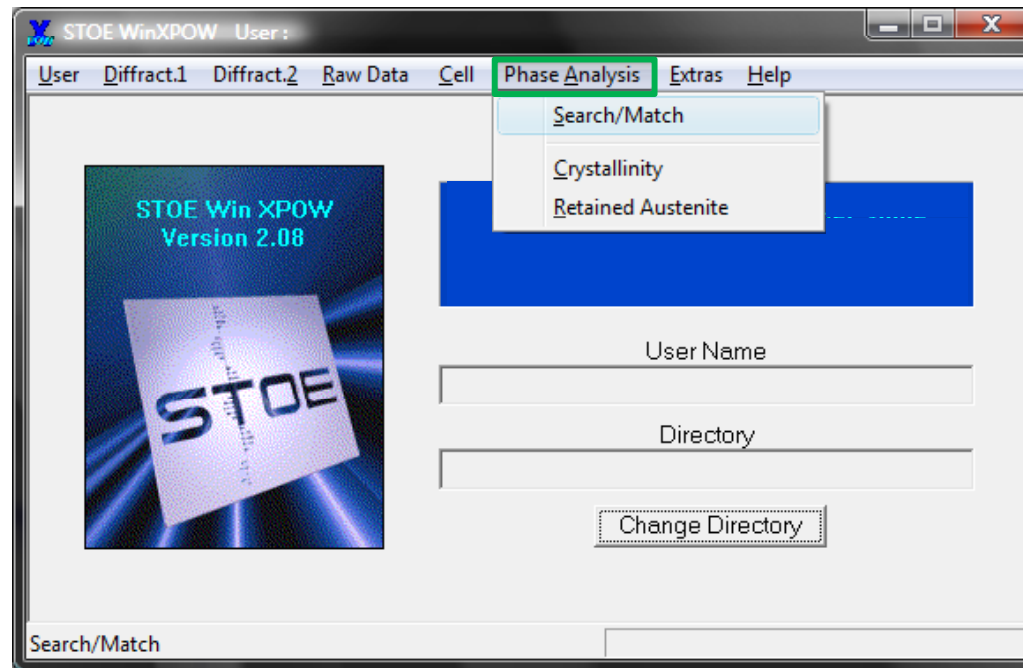


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Phase Analysis → Search Match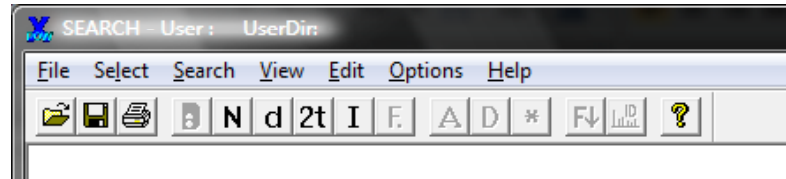

- Select → Elements

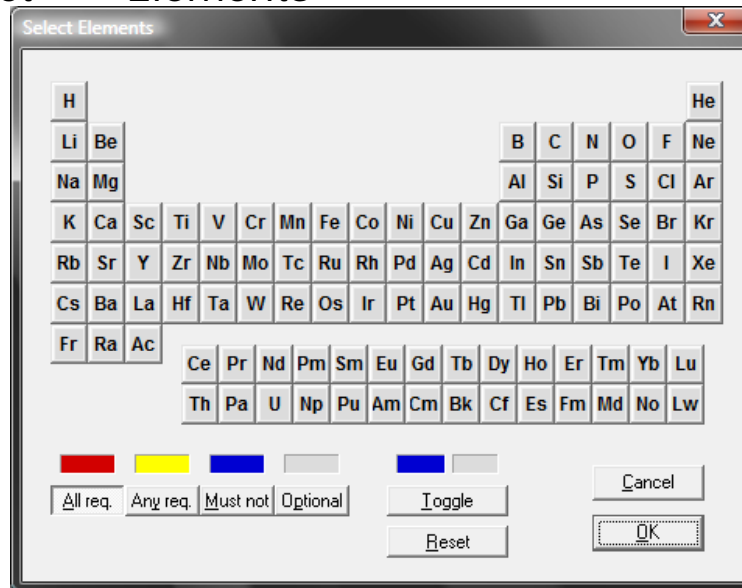
1. press **Toggle**


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Phase Analysis → Search Match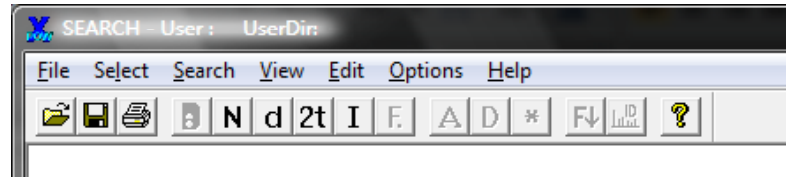

- Select → Elements

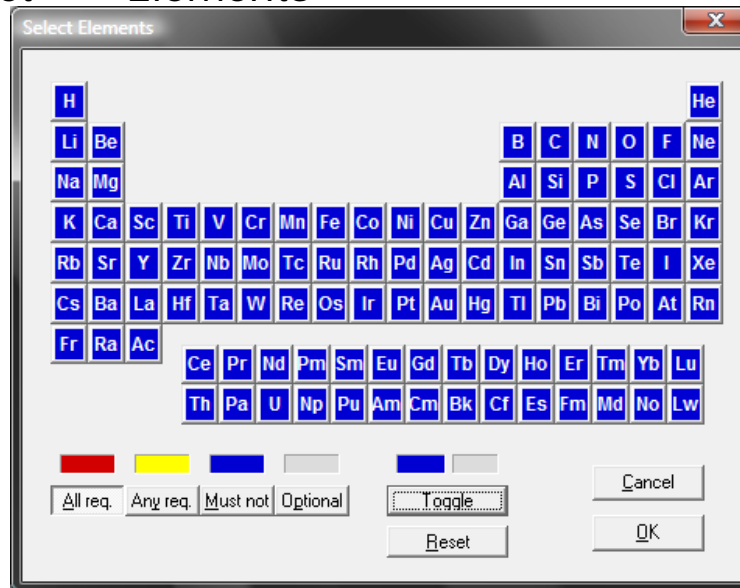
1. press **Toggle**
2. select **All required elements**


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Phase Analysis → Search Match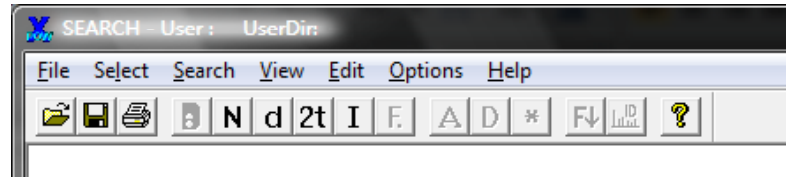

- Select → Elements

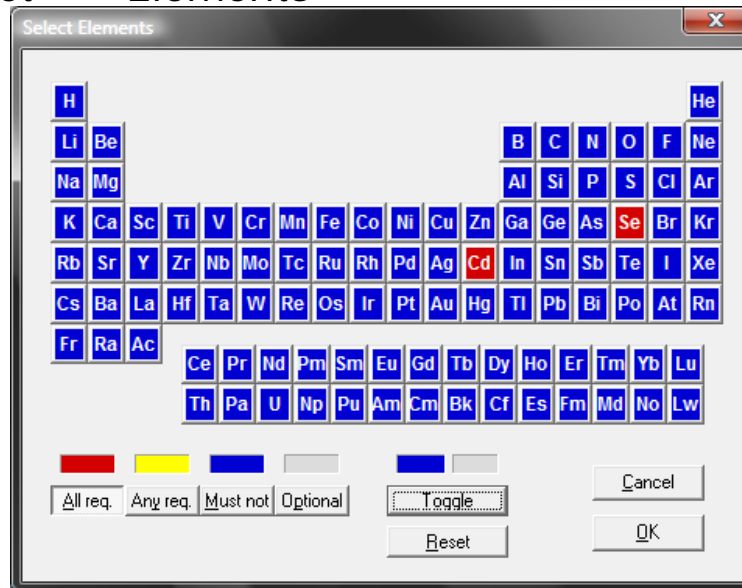
1. press **Toggle**
2. select **All required elements**
3. press **OK**


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Phase Analysis → Search Match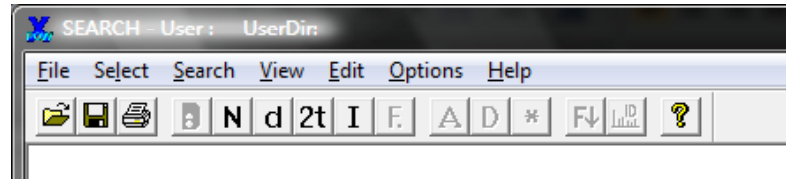

- Select → Elements

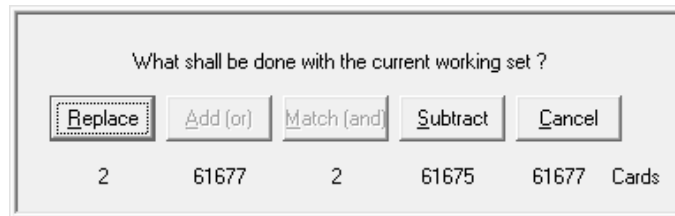1. press **Toggle**
2. select **All required elements**
3. press **OK**
4. press **Replace**

AK Eychmüller GroupSeminar


Brief introduction to the WinXPow for XRD.

- Phase Analysis → Search Match

view Names

File	Select	Search	View	Edit	Options	Help
						
1 [8- 459] Cd Se / Cadmium Selenide / Cadmoselite, syn						
2 [19- 191] Cd Se / Cadmium Selenide						


view d-values

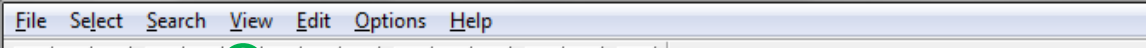File	Select	Search	View	Edit	Options	Help
						
1 [8- 459] 3.720(100) 3.510(70) 3.290(75) 2.554(35) 2.151(85)						
2 [19- 191] 3.510(100) 2.149(55) 1.833(25) 1.519(2) 1.394(4)						

view 2theta´s

File	Select	Search	View	Edit	Options	Help
						
1 [8- 459] 23.901(100) 25.353(70) 27.081(75) 35.108(35) 41.969(85)						
2 [19- 191] 25.353(100) 42.009(55) 49.698(25) 60.942(2) 67.085(4)						


view intensities

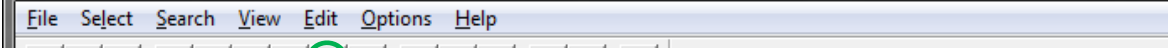File	Select	Search	View	Edit	Options	Help
						
1 [8- 459] 23.901(100) 41.969(85) 27.081(75) 25.353(70) 45.791(70)						
2 [19- 191] 25.353(100) 42.009(55) 49.698(25) 67.085(4) 76.779(4)						


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Raw Data → Graphics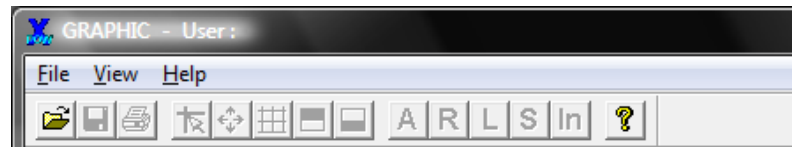

- File → Open → cw_Alloy-X1opt.raw

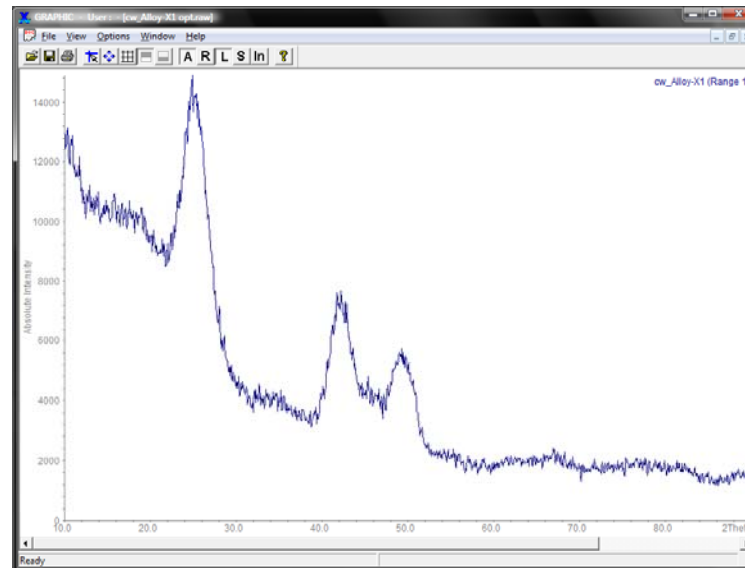AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Right Mouse Button:

Add Raw Data	to overlay different *.raw-files
Add Peak Data	to overlay a theoretical pattern or *.pks-file
Add ICDD Data	to overlay a theoretical pattern from the ICDD Data
Split	split the screen
Scaling Axes	scales the axis
Edit Curves ...	option to delete overlayed data
Insert Note ...	insert a note


AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Right Mouse Button → Add ICDD Data


type in the number with space or komma

→ OK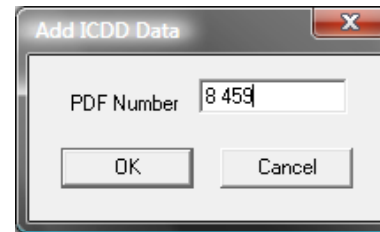

Right number? Correct substance?

→ Ja

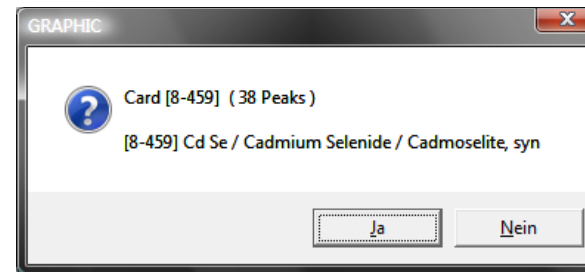AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Right Mouse Button → Add ICDD Data

Right number! Correct substance!

AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Right Mouse Button → Add ICDD Data

Right number! Correct substance!

AK Eychmüller GroupSeminar

Brief introduction to the WinXPow for XRD.

- Shut down WinXPow:

first:

close every working window (e.g. GRAPHICS, SEARCH/MATCH)

last:

close MAIN MENU